


Instructions for TEC controller

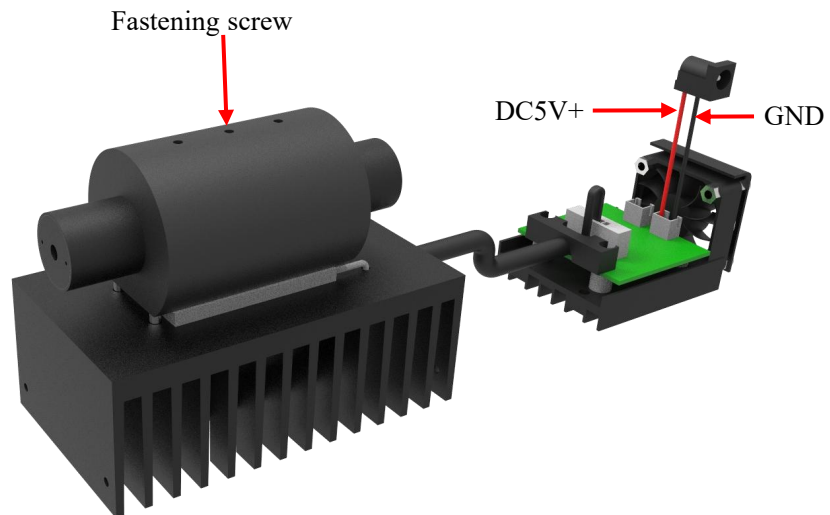

1. The red line on the power supply of the TEC controller is “+”, please connects it with 5VDC. The black line is “-”, please connect it with GND.
2. Loosen the fastening screw and insert the laser to the TEC controller, then tighten the fastening screw. After that take off the label pasted at the aperture of the laser module.
3. Connect the adapter with the earphone socket on the laser, then connect the adapter to external power supply and the laser emits. Waiting for the output power stable before use.
4. Closing the laser system: Turn off the mains power of the power supply.
5. To prevent optic path from dust, you should replace aperture label.